STANDAR OPERASIONAL PROSEDUR (SOP)
PENGUNAAN ALAT LABORATORIUM
(STANDARD OPERATING PROCEDURE (SOP)
FOR THE USE OF LABORATORY EQUIPMENT)

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The author expresses gratitude to Almighty God, by whose grace and blessings, the Standard Operating Procedure (SOP) for the Use of Chemistry Laboratory Equipment at FMIPA UHO has been successfully completed. The preparation of this SOP is intended to meet the need for standard operating procedures as an effort to improve the quality and performance of the laboratory in the future. This SOP serves as a guideline for students, lecturers, and staff in carrying out their daily duties. The author welcomes criticism and suggestions for the future perfection of this SOP in order to establish good administrative order. Thus, this SOP is prepared with the hope that it can assist in improving the management of the work environment at the Chemistry Laboratory of FMIPA UHO

Kendari,

Kepala Laboratorium Kimia FMIPA UHO

Head of Chemistry Laboratory FMIPA UHO

Dr. Sitti Hadijah Sabarawati, S.Si., M.Si.

NIP.19710514 200003 2 001

DAFTAR ISI

(TABLE OF CONTENTS)

Judul (Title)	Halaman (Page)
HALAMAN JUDUL (<i>TITLE PAGE</i>)	-
KATA PENGANTAR (<i>FOREWORD</i>)	1
DAFTAR ISI (<i>TABLE OF CONTENTS</i>)	2
I. Definisi (<i>Definition</i>)	3
II. Tujuan (<i>Objective</i>)	3
III. Ruang Lingkup (<i>Scope</i>)	3
IV. Definisi Istilah (<i>Definition of Terms</i>)	3
V. Prosedur Operasi Standar (SOP) Penggunaan Alat Laboratorium <i>Standard Operating Procedure (SOP) for the Use of Laboratory Equipment</i>	4

STANDARD OPERATING PROCEDURE (SOP) CHEMISTRY LABORATORY FMIPA UHO)

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The Chemistry Laboratory of FMIPA UHO is an academic facility of the Faculty of Mathematics and Natural Sciences UHO that provides services in the form of facility usage, equipment, consumables, and expertise for the purposes of practicum, research/community service, and services/analysis.

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The Standard Operating Procedure aims to provide guidelines for the process of utilizing the laboratory for practicum services, research, and services/analysis by users.

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The Chemistry Laboratory of FMIPA UHO provides services for students of FMIPA and other faculties within UHO, as well as the general public.

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Head of the laboratory is a lecturer from the Department of Chemistry who is given the additional task of leading and coordinating all activities taking place in the Chemistry Laboratory.

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Laboratory technician (Laboran) is a technician assigned to assist the Head of the Laboratory in organizing all activities in the Laboratory

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Practicum assistant is a student selected and appointed by the laboratory to assist the supervising lecturer in overseeing the implementation of the practicum conducted by the Practicum Students.

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Practicum student (Praktikan) is a student conducting practical work in the Analytical Chemistry Laboratory during the respective semester.

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Service user refers to students, lecturers, and external parties who utilize the Laboratory.

V. PROSEDUR OPERASI STANDAR (SOP) SOP PENGGUNAAN ALAT LABORATORIUM *(SOP FOR THE USE OF LABORATORY EQUIPMENT)*

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The oven is only used for drying practicum/research materials/substances that are non-corrosive.

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